



# **NMOS** multichannel detector heads

C5964 series

# UV to near infrared range (200 to 1000 nm), For multichannel spectrophotometry

The C5964 series is a family of multichannel detectors developed for spectrophotometry in the UV to near infrared range (up to 1000 nm). The C5964 series device incorporates a thermoelectrically-cooled NMOS linear image sensor (S5930/S5931/ S8382/S8383 series), low noise driver/amplifier circuit and highly stable temperature control circuit. It also operates from simple external signal inputs. The image sensor is cooled to a preset temperature (Ts=0 °C) as soon as the power is turned on. Should the cooler fail and the device overheat, the built-in protection circuit automatically shut off the power. The housing is designed for compactness, yet offers good heart dissipation. Furthermore, mounting holes are provided on the front panel of the housing, permitting easy connection to a monochromator or other instruments

#### Features

- Designed for use with a thermoelectrically-cooled NMOS linear image sensors (S5930 series, etc.)
- Built-in driver/amplifier and temperature control circuits
- Highly stable temperature control ensures a constant cooling temperature of Ts=0  $\pm$  0.05 °C (at Ta=10 to 30 °C)
- Operates from simple signal inputs
- High sensitivity and wide dynamic range

# Applications

- Multichannel spectrophotometry
- Spectrophotometer, colorimeter
- Optical spectrum analyzer
- **Time-resolved photometry**

#### Selection guide

The C5964 series consists of the following models depending on the NMOS linear image sensor used.

NMOS	NMOS linear image sensor					
multichannel	Type no	Type no Number of pivols		Effective image area	Remark	
detector head	Type no.	Number of pixels	[μm (H) × μm (V)]	[mm (H) × mm (V)]	Remark	
C5964-0800	S5930-256S	256	50 × 2500	12.8 × 2.5		
C5964-0900	S5930-512S	512	30 × 2300	25.6 × 2.5	Standard type	
C5964-0910	S5931-512S	512	25 × 2500	12.8 × 2.5		
C5964-1010	S5931-1024S	1024	25 × 2500	25.6 × 2.5		
C5964-0901	S8382-512S	512	50 × 2500	25.6 × 2.5	IR-enhanced type	
C5964-1011	S8383-1024S	1024	25 × 2500	25.6 × 2.5	in-eminanced type	

### **■** Absolute maximum ratings

Parameter	Symbol	Value	Unit
Supply voltage (for digital circuitry)	+VD	-0.5 to +7	V
Supply voltage (for analog circuitry)	±VA	±18	V
Digital input voltage	-	VD	V
Operating temperature	Topr	+10 to +30	°C
Storage temperature	Tstg	0 to +50	°C

Note: Exceeding the absolute maximum ratings even momentarily may cause a drop in product quality. Always be sure to use the product within the absolute maximum ratings.

# **■** Electrical characteristics (Ta=25 °C, VD=+5 V, ±VA=±15 V, unless otherwise noted)

Parameter			Symbol	Min.	Тур.	Max.	Unit
Digital input		High level	VIH	+2.0	-	+VD	V
		Low level	VIL	-0.5	-	+0.8	V
Master clock (CLK) pulse frequency		fclk	-	-	375	kHz	
Video signal readout frequency			fv	-	-	fclk/6	Hz
Master start (Start) pulse width			tst	1/fclk	-	-	S
Didital Olitolit		High level (Io=-6 mA)	VIH	+2.0	-	-	V
		Low level (Io=+6 mA)	VIL	-	-	+0.8	V
Power supply conditions  Rated voltage  Current consumption	Digital	+VD	+4.75	+5.0	+5.25	V	
	Rateu voitage	Analog	±VΑ	±14.5	±15.0	±15.5	V
	Current consumption	+5 Vdc*1	-	-	-	+2.0	Α
		+15 Vdc	-	-	-	+100	mA
		-15 Vdc	-	-	-	-100	mA

<sup>\*1:</sup> Including the current consumption of the Peltier element incorporated in the NMOS linear image sensors (S5930 series, etc.)

# **■** Electrical and optical characteristics (Ta=25 °C, Tchip=0 °C, VD=+5 V, ±VA=±15 V, unless otherwise noted)

Parameter	Built in sensor	Symbol	Min.	Тур.	Max.	Unit
Spectral response range		λ	-	200 to 1000	-	nm
Peak sensitivity wavelength	S5930/S5931 series	\n	-	600	-	nm
reak sensitivity wavelength	S8382/S8383 series	λр	-	750	-	
	S5930 series		-	50	-	pC
Caturation output charge	S5931 series	Ocat	-	25	-	
Saturation output charge	S8382 series	Qsat	-	50	-	
	S8383 series		-	25	-	
	S5930 series		-	0.009	0.03	- pA/pixel
Dark current	S5931 series	ID	-	0.004	0.01	
Dark Current	S8382 series		-	0.009	0.03	
	S8383 series		-	0.004	0.01	
Photoresponse nonuniformity*2		PRNU	-	-	±3	%
	S5930 series		-	0.2	-	- V/pC
Conversion gain	S5931 series	Co	-	0.4	-	
Conversion gain	S8382 series	Gc	-	0.2	-	
	S8383 series		-	0.4	-	

 $<sup>^{\</sup>star}2:$  Use for  $^{\backprime}\!A''$  light source. 50% of saturation, excluding first and last pixels.




### **Temperature controller specifications (Ta=25 °C, VD=+5 V, ±VA=±15 V, unless otherwise noted)**


Parameter*3	Symbol	Min.	Тур.	Max.	Unit
Cooling temperature	Ts	-1	0	+1	°C
Temperature control range	ΔTs	-0.05	-	+0.05	°C
Power dissipation of Peltier element	Рр	-	-	7	W
Cool down time to preset temperature	to	-	-	5	min
Setting temperature for overheart protection	То	+40	-	-	°C

<sup>\*3:</sup> Other functions include error display, automatic power off, and detection of electrical opens and shorts by the thermosensor.

# Spectral response (typical example)

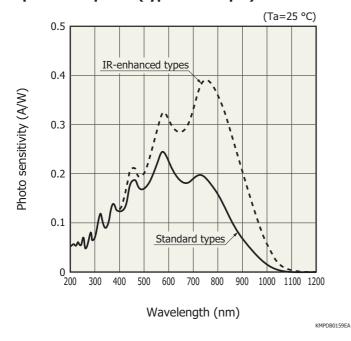


# **Equivalent circuit (NMOS linear image sensor S5930/S5931/S8382/S8383 series)**

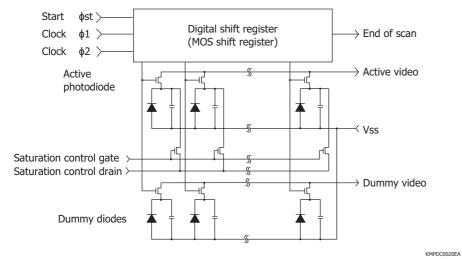

IAMATSU


### Block diagram (C5964 series)

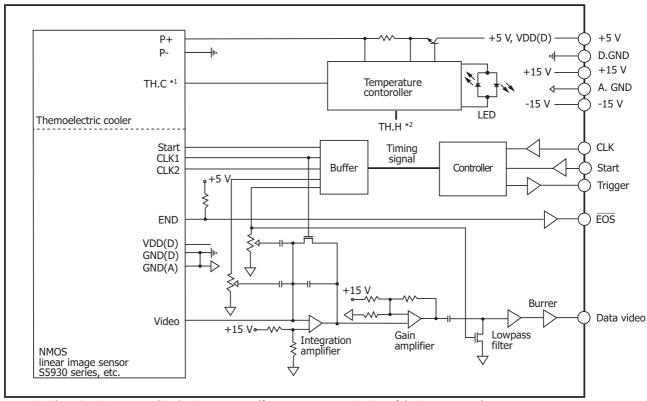


\*1: Thermistor incorporated in the image sensor (for temperature monitoring of the image sensor) \*2: Thermistor mounted on the heatsink fins (for temperature monitoring of the heat radiating side)

KACCC0067EA

## **Timing chart**

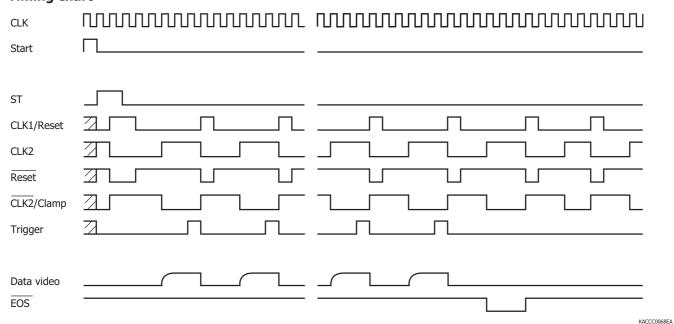


# Dimensional outline (unit: mm)

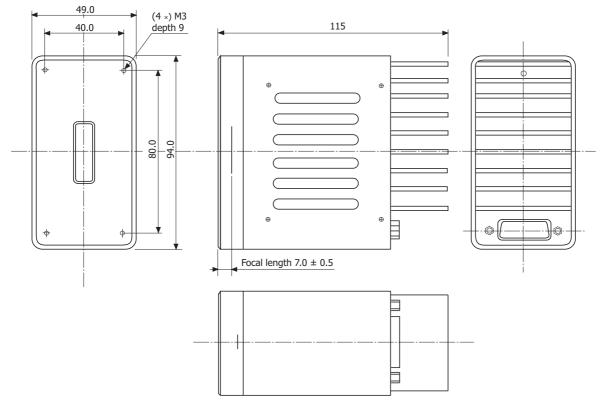

KACCA0053EA

#### **Pin connections**

15-pin D-sub connector (socket type)

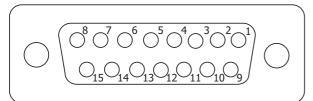

KACCC0069EA

Pin no.	Symbol	Description
1	NC	No connection
2	Data video	Analog video output signal. Positive polarity
3	VA1+(+15 V)	Power supply for analog circuitry
4	Va1-(-15 V)	Power supply for analog circuitry
5	VD(+5 V, P+)	Power supply for digital circuitry. For the thermoelectric cooler in the NMOS linear image sensor
6	Start	Digital input signal to initialize the circuit HCMOS compatible. Positive logic The start pulse interval determines the signal storage time of the sensor.
7	CLK	Digital input signal to specify the circuit operation HCMOS compatible. Operates at the rising edge
8	EOS	Digital output signal to indicate the end of scan of the NMOS linear image sensor HCMOS compatible. Negative logic
9	A.GND	Analog ground
10	A.GND	Analog ground
11	NC	No connection
12	D.GND(P-)	Digital ground. Power supply return of the thermoelectric cooler mounted in the NMOS linear image sensor
13	D.GND	Digital ground
14	D.GND	Digital ground
15	Trigger	Digital output signal for A/D conversion HCMOS compatible. Positive logic

# Precautions

The C5964 series is a precision device, so use sufficient caution when handling it.


- · Never disassemble or modify the device as this may cause an operating failure.
- · Protect the device from shocks such as drops or impacts.
- · Avoid storing the device in high temperature and high humidity locations for long periods of time.
- $\cdot \text{ Never block the air vents provided on the top and side panels of this unit during operation. Blocking the air vents may cause overheating.}$
- · When connecting to other equipment, check that the electrical specifications are matched.
- · Never exceed the maximum ratings during operation.

Observe the following precautions to obtain the fullest performance of device.

- · Provide a proper shield to protect the device against external electromagnetic effects. Use of a shield cable is recommended.
- $\cdot$  Use a power supply with minimum ripple and noise.
- · Pay attention to prevent extraneous light from entering the device for accurate measurements.



#### Connection example

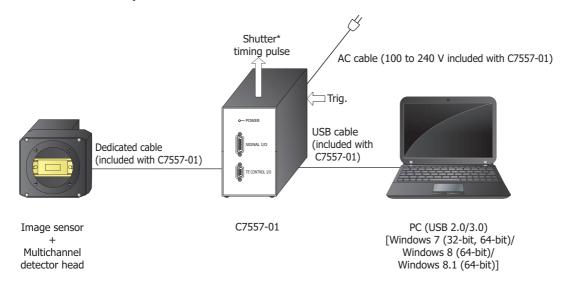

<sup>\*</sup> Shutter, etc. are not available.

KACCC0402FD

### **Precaution for use**

When operating the C5964 series with the C7557-01, always be sure to attach the MOS adapter (supplied) to the C7557-01 main unit. If the the C5964 series is connected and the power is turned on without attaching the adapter, the power supply in the C7557-01 main unit may be damaged.

#### Related information

www.hamamatsu.com/sp/ssd/doc\_en.html

- Precautions
- Disclaimer

Information described in this material is current as of October 2016.

Product specifications are subject to change without prior notice due to improvements or other reasons. This document has been carefully prepared and the information contained is believed to be accurate. In rare cases, however, there may be inaccuracies such as text errors. Before using these products, always contact us for the delivery specification sheet to check the latest specifications.

The product warranty is valid for one year after delivery and is limited to product repair or replacement for defects discovered and reported to us within that one year period. However, even if within the warranty period we accept absolutely no liability for any loss caused by natural disasters or improper product use. Copying or reprinting the contents described in this material in whole or in part is prohibited without our prior permission.

www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Solid State Division

1126-1 Ichino-cho, Higashi-ku, Hamamatsu City, 435-8558 Japan, Telephone: (81) 53-434-3311, Fax: (81) 53-434-5184

1126-1 ICnino-Cho, Higasni-Ku, Hamamatsu City, 435-8558 Japan, Telephone: (81) 53-434-3511, FaX: (81) 53-434-5184
U.S.A.: Hamamatsu Corporation: 360 Foothill Road, Bridgewater, N.J. 08807, U.S.A., Telephone: (1) 908-231-0960, Fax: (1) 908-231-1218
Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerst: 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49) 8152-375-0, Fax: (49) 8152-265-8
France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: 33-(1) 69 53 71 00, Fax: 33-(1) 69 53 71 00
United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road, Welwyn Garden City, Hertfordshire AL7 1BW, United Kingdom, Telephone: (44) 1707-294888, Fax: (44) 1707-32577
North Europe: Hamamatsu Photonics Norden AB: Torshamnsgatan 35 16440 Kista, Sweden, Telephone: (46) 8-509-031-01
Italy: Hamamatsu Photonics Italia S.r.l.: Strada della Moia, 1 int. 6, 20020 Arese (Milano), Italy, Telephone: (39) 02-93581733, Fax: (39) 02-93581741
China: Hamamatsu Photonics (China) Co., Ltd.: B1201, Jiaming Center, No.27 Dongsanhuan Beilu, Chaoyang District, Beijing 100020, China, Telephone: (86) 10-6586-6006, Fax: (86) 10-6586-2866