

NEW

Scientific CMOS board level camera C11440-62U



The C11440-62U is a compact and highly sensitive board level camera with a scientific CMOS image sensor. The camera has over 80 % quantum efficiency at 600 nm. The camera runs at 30 frames/s at full resolution and it can reach up to 25 655 frames/s (512 × 8 ROI). The C11440-62U can acquire a high-quality image at longer exposure time by having an element cooling function built-in. The C11440-62U has low readout noise of 2.1 electrons (rms) at 30 frames/s. The compact (92 mm × 92 mm) and light weight design is ideal for integration into various measurement and analytical instruments. In addition to that the camera employs a USB 3.0 digital interface for convenience and compatibility.


FEATURES

- Sensor temperature stabilization
- Resolution: 4.0 megapixels
- Readout noise: 2.1 electrons (rms)
: 1.5 electrons (median)
- Frame rate: 30 frames/s (Full resolution)
- Quantum efficiency: 82 % (Peak QE)


APPLICATIONS

- Super-resolution microscope
- Real time confocal microscope
- Live cell imaging
- Fluorescence time lapse imaging
- DNA sequence
- DNA microarray
- X-ray I.I., X-ray scintillator readout

SPECTRAL RESPONSE

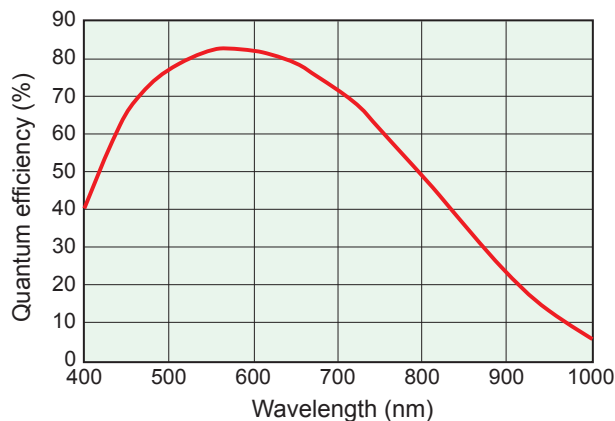

SYSTEM CONFIGURATION

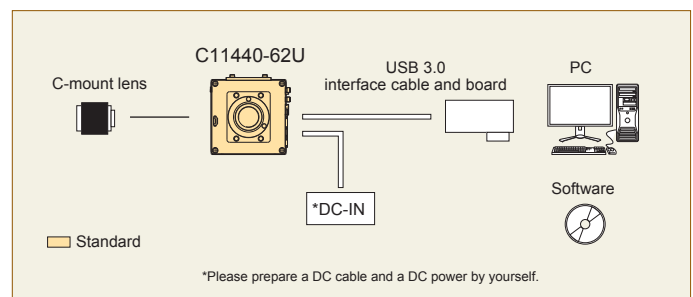

SPECIFICATIONS

Type number	C11440-62U
Imaging device	Scientific CMOS image sensor
Effective number of pixels	2048 (H) × 2048 (V)
Cell size	6.5 μm (H) × 6.5 μm (V)
Effective area	13.312 mm (H) × 13.312 mm (V)
Full well capacity typ.	30 000 electrons
Dynamic range typ.*1	20 000 : 1
Frame rate	30 frames/s (Full resolution)
	25 655 frames/s (512×8 ROI)
Readout noise typ.*2	2.1 electrons (rms), 1.5 electrons (median)
Digital output	16 bit
Exposure time *3	1 ms to 30 s (Internal trigger mode at full resolution)
External trigger input mode	Edge trigger, Level trigger, Synchronous readout trigger, Start trigger, Global reset edge trigger, Global reset level trigger
Trigger output	Programmable timing outputs, Low output, High output, Trigger ready output, Global exposure timing output
Binning *4	2×2, 4×4
Sub-array	Yes
Interface	USB 3.0
Lens mount	C-mount
Connector	USB 3.0 microB type
Power supply	DC+12 V
Power consumption	Approx.15 VA
Ambient operating temperature	0 °C to +40 °C
Ambient operating humidity	30 % to 80 % (with no condensation)
Ambient storage temperature	-10 °C to +50 °C
Ambient storage humidity	90 % max. (with no condensation)


*1: Calculated from the ratio of the full well capacity and the average readout noise.

*2: Value of the sensor temperature 27 °C at 33 ms exposure.

*3: When exposure time is longer than 100 ms, white spots would be visible and image quality would be degraded.

*4: Digital binning processing in the camera

[Sensor temperature stabilization function details]

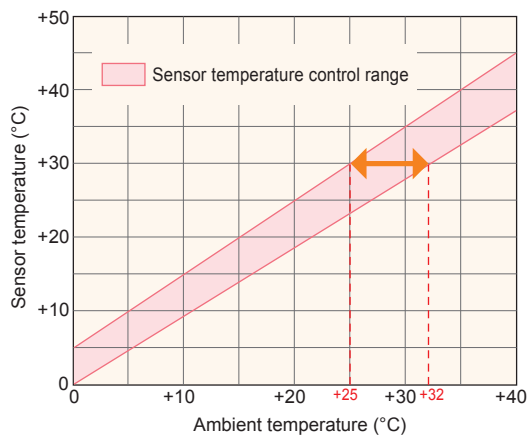

It is a range where the sensor temperature can be kept constant even if environmental temperature fluctuates.


Reference example (shown the orange arrow in figure):

The sensor temperature can be kept at +30 °C within the range of ambient temperature +25 °C to +32 °C.

- Product and software package names noted in this documentation are trademarks or registered trademarks of their respective manufacturers.
- Subject to local technical requirements and regulations, availability of products included in this promotional material may vary. Please consult your local sales representative.
- Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications and external appearance are subject to change without notice.

© 2018 Hamamatsu Photonics K.K.

DIMENSIONAL OUTLINES (Unit : mm)

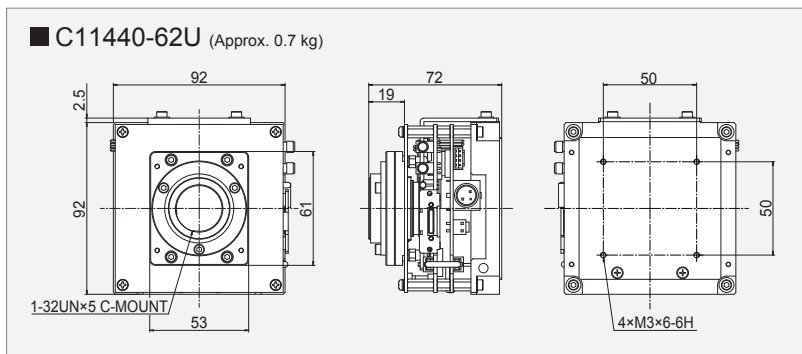

OPTIONS

- External trigger cable SMA-BNC 5 m: A12106-05
- External trigger cable SMA-SMA 5 m: A12107-05

HAMAMATSU PHOTONICS K.K. www.hamamatsu.com

HAMAMATSU PHOTONICS K.K., Systems Division

812 Joko-cho, Higashi-ku, Hamamatsu City, 431-3196, Japan, Telephone: (81)53-431-0124, Fax: (81)53-433-8031, E-mail: export@sys.hpk.co.jp

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, Bridgewater, NJ 08807, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: usa@hamamatsu.com

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-265-8 E-mail: info@hamamatsu.de

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: infos@hamamatsu.fr

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road, Welwyn Garden City, Hertfordshire AL7 1BW, UK, Telephone: (44)1707-294888, Fax: (44)1707-325777 E-mail: info@hamamatsu.co.uk

North Europe: Hamamatsu Photonics Norden AB: Torshamnsgatan 35 16440 Kista, Sweden, Telephone: (46)8-509 031 00, Fax: (46)8-509 031 01 E-mail: info@hamamatsu.se

Italy: Hamamatsu Photonics Italia S.r.l.: Strada della Moia, 1 int. 6, 20020 Arese (Milano), Italy, Telephone: (39)02-93 58 17 33, Fax: (39)02-93 58 17 41 E-mail: info@hamamatsu.it

China: Hamamatsu Photonics (China) Co., Ltd.: 1201 Tower B, Jiaming Center, 27 Dongsanhuan Beilu, Chaoyang District, 100020 Beijing, China, Telephone: (86)10-6586-6006, Fax: (86)10-6586-2866 E-mail: hpc@hamamatsu.com.cn

Taiwan: Hamamatsu Photonics Taiwan Co., Ltd.: 8F-3, No.158, Section2, Gongdao 5th Road, East District, Hsinchu, 300, Taiwan R.O.C. Telephone: (886)03-659-0080, Fax: (886)03-659-0081 E-mail: info@hamamatsu.com.tw

Cat. No. SCAS0115E03
JUN/2018 HPK
Created in Japan